


No. 4

Schlieren Imaging of Shock Wave from Trigger Air Spark Gap

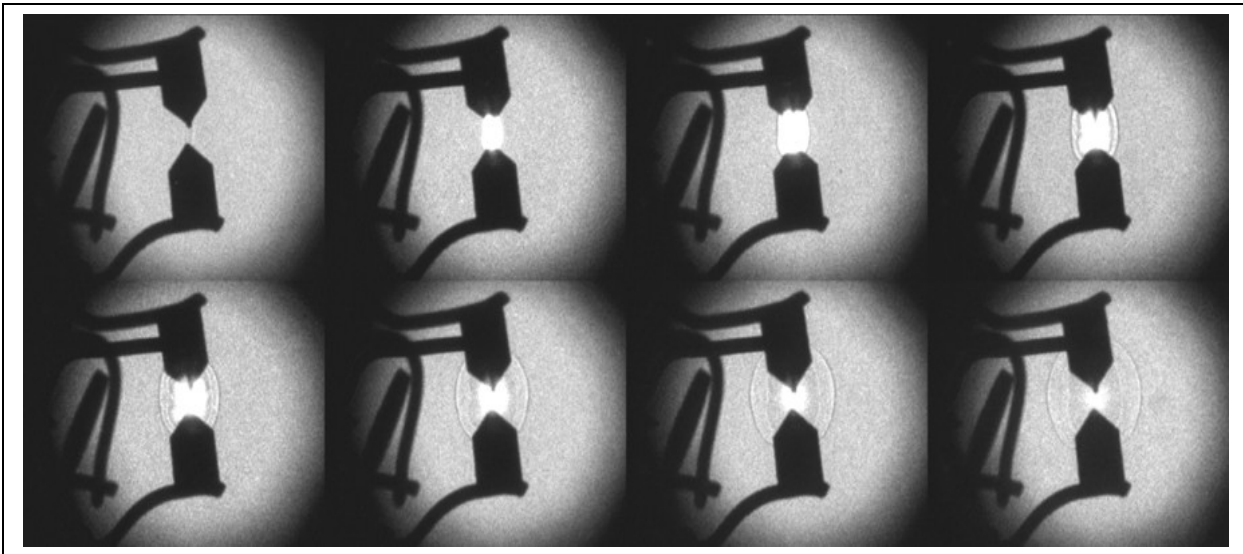

IMAGING PARAMETERS

SIM8 Camera running at 5 Million fps, 35 ns
Exposures with Pseudo Schlieren Optical system. The
field of view is about 10cm x 8cm.

OVERVIEW OF EXPERIMENT

The Air Spark Gap with 0.75 uF Capacitor and 1.5KV
Charge Voltage across the 2 mm air gap.
Two of the SIM's output triggers were used to synchroniz
the triggering of the spark gap and to pre-trigger the flash
lamp used for the Schlieren light source.

EQUIPMENT PARAMETERS

A 1.5 msec duration xenon flash lamp directly coupled to a 0.125 inch diameter fibre light guide which
was the point source. A 200 mm f2.5 lens was used to collimate the point source into a parallel beam,
The Objective lens, a Nikon 85 mm f1.8 with extension tubes, was used to both focus the event into
the camera at about 1:1 magnification. With the lens position away from the camera using the
extension tube, the parallel light focuses at the infinity focus point of the objective lens, which then re-
diverges to produce the background light on the image plane to produce the Schlieren imaging effect.

UK (Head Office / Factory)

6 Harvington Park, Pitstone Green
Business Park, Pitstone.
LU7 9GX England

+44 (0) 1442 827728

USA

Specialised Imaging Inc.
40935 County Center Dr. Suite D
Temecula, CA 92591, USA

+1 951-296-6406

GERMANY

Hauptstr. 10,
82275 Emmering
Germany

+49 8141 666 89 50



FM 87429 ISO9001:2015

